

ScienceCube



# Wireless Salinity (WL119S) User Guide



Rev. WL119S-12-2023

**This product is to be used for educational purposes only. It is not appropriate for industrial, medical, research, or commercial applications.**

 **KOREADIGITAL**

*The Science Cube wireless Salinity sensor can measure the conductivity of solution.*

**The wireless salinity sensors** can measure the amount of salt dissolved in water. Salinity can be measured in a variety of environments, from plain water to seawater. The salinity of sea water is about 35ppt.

You can measure by remotely connecting to a smart device or PC wirelessly or wired.

## Suggested experiments

Using a wireless salinity sensor, you can perform experiments such as:

- Measurement of salinity of seawater
- Seasonal salinity measurement
- Salinity measurement in aquarium
- Comparison of salt water concentration percentages

## Composition

*The ScienceCube wireless conductivity sensor consists of the following.*

- Wireless salinity sensor(WL119S)
- USB-A/C cable
- Booklet

## Feature

- Up to four Science Cube wireless sensors can be connected to a PC or smart device at the same time.
- It supports dual-mode Bluetooth, allowing you to connect not only smart devices

but also desktop and laptop PCs to conduct experiments using the **Science#** application.

- It can be connected to a PC through a USB port and experiments can be performed using the **Science#** program.



## Function of wireless sensor

### Structure




- ① USB port : Connect the sensor to a PC and use it for experiments or charging.
- ② OLED Display : Displays measured sensor values, sensor type, sensor ID, and remaining battery level.
- ③ Power/Function Button : It has functions such as power ON/OFF, measurement sensor change and calibration, etc.
- ④ Sensing part : A part immersed in the solution to be measured, where opposing

electrodes and a temperature probe are located.

### Power/Function Button

Status	Turn	Action	Description
When the power is off	Click once	■	A short press turns the sensor on.
	Long click	■■■■■■■■■■	A long press changes the mode and turns on the sensor.
When it's on	Double click	■■■	Zero point setting (For sensors with zero point function) is performed and <b>U0</b> is displayed above the device.
	Long click	■■■■■■■■■■	Turns off.


### Start screen

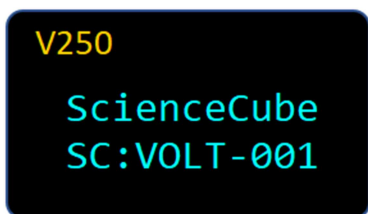

V250 : Displays the sensor's firmware version.

SC:XXXX-001 : When you search for a Bluetooth device, the device name will be displayed. (Sensor name and 3-digit serial number)

### Mode change

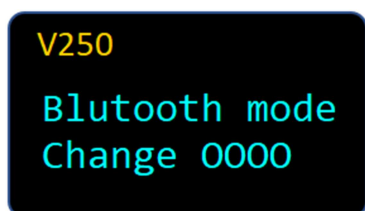

When you press and hold the power button and turn it on, the Bluetooth connection mode changes to **Mobile** or **PC** with the following message.

## Measurement screen



<p>① <b>Connection mode</b></p>	<p><b>Mobile</b> : Connecting Android or iOS.  <b>PC</b> : Connecting to Windows PC          ※ A long press changes the mode and turns on the sensor.</p>
<p>② <b>Sensor-ID</b></p>	<p>This is the sensor's unique number and is displayed along with the sensor name in the device name when connected via Bluetooth.</p>
<p>③ <b>Battery</b></p>	<p>Check the battery status, and when charging via USB, the display will change to charging.</p>
<p>④ <b>Value</b></p>	<p>1) Displays sensor measurement values and units in real time.          2) If <b>user calibration</b> is used, <b>U0</b> or <b>UC</b> will be displayed above the units.          3) For sensors with <b>multiple ranges</b>, the current range is displayed.          4) For <b>multiple sensors</b>, the values for each sensor type are displayed.</p>

## How it Works

A salinity sensor is a sensor that measures the salinity of water. Salinity is an indicator of the amount of salt dissolved in water. Salinity sensors calculate salinity by measuring electrical conductivity. In the case of seawater, salinity and electrical conductivity are proportional, so salinity can be easily calculated by measuring electrical conductivity.

ScienceCube salinity sensor measures the ability of a solution to conduct an electric current between two electrodes. In solution, the current flows by ion transport. Therefore, an increasing concentration of ions in the solution will result in higher conductivity values. The salinity probe is actually measuring conductance, defined as the reciprocal of resistance.

Conductivity,  $C$ , is found using the following formula:


$$C = G \cdot kc$$

where  $G$  is the conductance, and  $kc$  is the cell constant. The cell constant is determined

for a probe using the following formula:

$$kc = d / A$$

where  $d$  is the distance between the two electrodes, and  $A$  is the area of the electrode surface. The conductivity value is found by multiplying conductance and the cell constant. Since the ScienceCube® Salinity Probe also has a cell constant of  $10 \text{ cm}^{-1}$ .

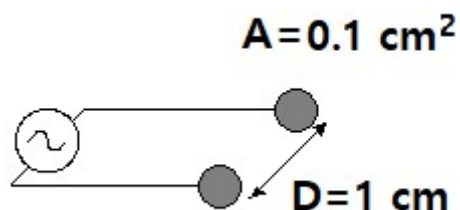

## Using the Sensor

- Rinse the tip of the probe with distilled water. **Optional** : Blot the inside of the electrode cell dry only if you are concerned about water droplets diluting or contaminating the sample to be tested.
- Insert the tip of the probe into the sample to be tested. **Important** : Be sure the electrode surfaces in the elongated cell are completely submerged in the liquid.
- While gently swirling the probe, wait for the reading on your sensor to stabilize. This should no more than 5 to 10 seconds.
- Rinse the end of the probe with distilled water before taking another measurement.
- If you are taking readings at temperatures below 15°C or above 30°C, allow more time for the temperature compensation to adjust and provide a stable conductivity reading.


**Important** : Do not place the electrode in viscous, organic liquids, such as heavy oils, glycerin(glycerol), or ethylens glycol. Do not place the probe in acetone or non-polar solvents, such as pentans or hexane.

## Calibration

All sensors are precisely calibrated using standard solutions during the manufacturing process before being shipped and are ready for immediate use. However, you can perform **[User Calibration]** for more accurate measurements.

**[User Calibration]** is reflected temporarily and then returns to the default when the sensor is turned on again.

1. Press the Power/Calibrate button twice in succession.
2. When the sensor shows 0.00 ppt.
3. **[User Calibration]** is complete and the word "U0" is displayed above the unit on the sensor screen.

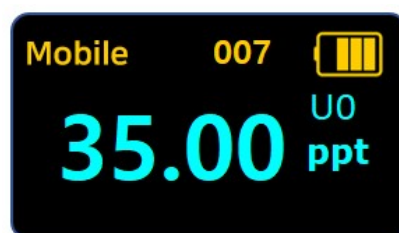

To cancel **[User Calibration]**, press the button twice in succession. U0 disappears and returns to default.

## Automatic Temperature Compensation

ScienceCube wireless salinity sensor is automatically temperature compensated between temperatures of 5 and 35°C. Note that the temperature of a solution is being read by a thermistor that extends into the space between the graphite electrodes. Readings are automatically referenced to a conductivity value at 25°C-therefore the Salinity Probe will give the same conductivity reading in a solution that is at 15°C as it would if the same solution were warmed to 25°C. This means you can calibrate your probe in the lab, and then use these stored calibrations to take readings in colder (or warmer) water in a lake or stream. If the probe was not temperature compensated, you would notice a change in the conductivity reading as temperature changed, even though the actual ion

concentration did not change.

## Using the Salinity sensor with Other ScienceCube Sensors

It is important to know what interaction occurs when you wire a wireless salinity sensor to a PC or smart device via a USB cable and measure with another wired sensor in the same solution (in the same aquarium or beaker, for example).

This situation arises because the Salinity Probe outputs a signal in the solution, and this signal can affect the reading of another probe.

**Note :** This does not apply if each device is connected wirelessly.

The sensors below cannot be used wired to the same PC or smart device as conductivity probes from the same solution:

- Dissolved Oxygen Probe
- pH or ORP Probe and Ion Selective Electrodes

## Storage and maintenance

- When you have finished using the Salinity Probe, simply rinse it off with distilled water and blot it dry using a paper towel or lab wipe. The probe can then be stored dry.
- If the probe cell surface is contaminated, soak it in water with a mild detergent for 15 minutes. Then soak it in a dilute acid solution (0.1 M hydrochloric acid or 0.5 M acetic acid works well) for another 15 minutes. Then rinse it well with distilled water. Important: Avoid scratching the inside electrode surfaces of the elongated cell.

**CAUTION: The electrode surface is treated with graphite. Do not touch or contaminate this area.**

## Specifications

Item	Description
Range	0 ~ 50 ppt
Resolution	0.01 ppt
Sampling Time	Max. 100Hz (0.01 sec.), (Typical 1Hz)
Condition	5 ~ 35°C (Automatic Compensation)
Wireless Connection	Bluetooth 5.0 or Classic 2.1
Wired Connection	USB-C
Battery	700mAh Li-Polymer rechargeable
Charging Time	within 2 hours
Operating Time	Approximately 6 hours after full charge (depending on usage conditions)

**CAUTION: Do not use the instrument beyond the measurement range or in conditions that exceed the short-term exposure limits. Prolonged exposure beyond the maximum permissible range can cause serious damage to the sensor.**

- ScienceCube® is a registered trademark of Korea Digital. Science# is a trademark of Korea Digital. All other trademarks are the property of their respective owners.
- The copyright of all products (hardware, software, content) related to Science Cube belongs to Korea Digital Co., Ltd.
- The contents of this manual are provided for informational purposes only, and product specifications and functions may be changed without prior notice to improve performance.
- This product is designed for science education. No warranty is provided and no liability is assumed for errors in industrial testing or manufacturing process controls, medical analysis or controls, or commercial design applications.

**Contact us**

**TEL : +82-2-2109-8839**

**FAX : +82-2-2109-8881**

**[www.sciencecube.com](http://www.sciencecube.com)**

**Korea Digital Co., Ltd.**

#804 Ace Twin Tower 273 Digital-ro Guro-gu Seoul 08381 Korea

[www.koreadigital.com](http://www.koreadigital.com)